

Product Information

Leupeptin Ready Made Solution

Non-freezing solution, 5 mg/mL

SAE0153

Product Description

Storage Temperature: -20 °C

Leupeptin, also known as *N*-acetyl-L-leucyl-L-leucyl-L-argininal, is a tripeptide produced by *Actinomycetes*/*Streptomyces*. Leupeptin is a reversible competitive inhibitor of cysteine proteases and serine proteases.¹ Leupeptin acts by covalent binding to, respectively:²

- Catalytic cysteines of cysteine proteases
- Catalytic series of serine proteases

Leupeptin was first isolated from microbial sources as a mixture of two very similar forms:³

- Acetyl-Leu-Leu-Arg-al
- Propionyl-Leu-Leu-Arg-al

While the propionyl leupeptin is active as an inhibitor, the acetyl form is more commonly used.

Leupeptin has been reported to inhibit calpain,⁴ cathepsin B,⁵ cathepsins H and L,⁶ and trypsin.⁷

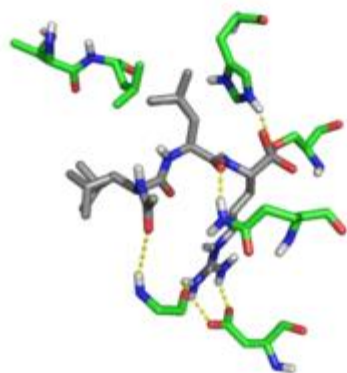


Figure 1. Crystal structure of Leupeptin (silver) in the Trypsin (green) binding pocket. Hydrogen bonds are shown as yellow dotted lines.^{8,9}

A typical working concentration range is 10-100 μ M. The activity of leupeptins and related analogs has been studied.¹⁰

HPLC analysis of leupeptin gives multiple peaks because of the formation of tautomeric isomers in aqueous solution.¹¹ The purity is determined using three main peaks. The majority of 'contaminating' peptide is racemized leupeptin.

Reagent

This Leupeptin Ready Made Solution is supplied as a proprietary non-freezing formulation, at 10 mM concentration (~5 mg/mL).

Precautions and Disclaimer

For R&D use only. Not for drug, household, or other uses. Please consult the Safety Data Sheet for information regarding hazards and safe handling practices.

Storage/Stability

The solution is shipped on cooler packs ("wet ice").

Storage at -20 °C is recommended. The product, as supplied, is stable for two years. For short time periods, the product can be stored at 2-8 °C.

Usage

The stock solution may be diluted 1:100 to obtain a 100 μ M working concentration. The effective concentration of leupeptin is in the range of 10 μ M-100 μ M.

Because of its aldehyde group, leupeptin may act as a reducing agent, and thus may interfere in protein determination assays, such as the Lowry assay and, to a lesser extent, the Bradford assay.

References

1. Umezawa, H., *Methods Enzymol.*, **45**, 678-695 (1976).
2. Kędzior, M. *et al.*, *Med. Microbiol. Immunol.*, **205(4)**, 275-279 (2016).
3. Aoyagi, T. *et al.*, *J. Antibiot. (Tokyo)*, **22(6)**, 283-286 (1969).
4. Zimmerman, U.-J.P. and Schlaepfer, W.W., *Biochemistry*, **21(17)**, 3977-3983 (1982).
5. Knight, C.G., *Biochem. J.*, **189(3)**, 447-453 (1980).
6. Zollner, H., *Handbook of Enzyme Inhibitors*, 2nd ed., Part B. VCH Press, pp. 821-822 (1993).
7. Kuramochi, H. *et al.*, *J. Biochem.*, **86(5)**, 1403-1410 (1979).
8. "1JRS: Hemiacetal Complex Between Leupeptin and Trypsin". RCSB PDB (Protein Data Bank), Released 14 October 1996 (<https://www.rcsb.org/structure/1JRS>).
9. Kurinov, I.V., and Harrison, R.W., *Protein Sci.*, **5(4)**, 752-758 (1996).
10. Maeda, K. *et al.*, *J. Antibiotics*, **24(6)**, 402-404 (1971).
11. Saino, T. *et al.*, *Chem. Pharm. Bull.*, **30(7)**, 2319-2325 (1982).

Notice

We provide information and advice to our customers on application technologies and regulatory matters to the best of our knowledge and ability, but without obligation or liability. Existing laws and regulations are to be observed in all cases by our customers. This also applies in respect to any rights of third parties. Our information and advice do not relieve our customers of their own responsibility for checking the suitability of our products for the envisaged purpose.

The information in this document is subject to change without notice and should not be construed as a commitment by the manufacturing or selling entity, or an affiliate. We assume no responsibility for any errors that may appear in this document.

Technical Assistance

Visit the tech service page at [SigmaAldrich.com/techservice](https://www.sigmaaldrich.com/techservice).

Standard Warranty

The applicable warranty for the products listed in this publication may be found at [SigmaAldrich.com/terms](https://www.sigmaaldrich.com/terms).

Contact Information

For the location of the office nearest you, go to [SigmaAldrich.com/offices](https://www.sigmaaldrich.com/offices).

The life science business of Merck operates as MilliporeSigma in the U.S. and Canada.

Merck and Sigma-Aldrich are trademarks of Merck KGaA, Darmstadt, Germany or its affiliates. All other trademarks are the property of their respective owners. Detailed information on trademarks is available via publicly accessible resources.

© 2020-2022 Merck KGaA, Darmstadt, Germany and/or its affiliates. All Rights Reserved.
SAE0153pis Rev 07/22 EO,CKV,DT,SM,GCY,MAM

The Merck logo, consisting of the word "MERCK" in a bold, red, sans-serif font.