

Product Information

Sephadex® G-25

Fine

G2580

Product Description

Sephadex® is a beaded gel filtration medium prepared by cross-linking dextran with epichlorohydrin under alkaline conditions.¹ General information and procedures for using gel filtration to separate proteins or to desalt protein solutions have been described.^{2,3}

This product can also be used for the separation of double-stranded DNA fragments. The exclusion limits for double-stranded DNA are as follows:

- G-25: 10 base pairs
- G-50: 20 base pairs
- G-100: 25 base pairs

DNA grade Sephadex® is available as part of our molecular biology product line, in the following Cat. Nos.:

- S5772 (G-25 Superfine)
- S5897 (G-50 Fine)
- S6022 (G-50 Medium)
- S6147 (G-100)

Cat. No. G2580 is an exact replacement for past Cat. Nos. 84942 and 27107-1.

Preparation Instructions

- This product should be placed in the usage buffer and allowed to swell for at least 3 hours at 20 °C or 1 hour at 90 °C.
- Once separation of the sample is complete, the gel should be washed with 2 column volumes of 0.2 M NaOH or a solution of non-ionic detergent, rinsed with water, and re-equilibrated with 2-3 column volumes of buffer.
- For storage, antimicrobial agents (such as 20% ethanol or 0.02% sodium azide) should be added to the suspension to prevent contamination.

- When necessary, the gel can be removed from the column and sterilized by autoclaving.
- The washed resin may then be resuspended in excess water or starting buffer, to pack the column bed.

Storage/Stability

Sephadex® does not melt and may be sterilized in its wet form at neutral pH by autoclaving for 30 minutes at 120 °C. This will not affect its chromatographic properties. However, if dry Sephadex® is heated to more than 120 °C, it will start to caramelize.

Sephadex® is stable in water, salt solutions, and organic and denaturing solvents. The pH stability is limited to low ionic strengths and short times when at the pH extremes of 2 and 13, particularly in the acid range. At low pH, partial hydrolysis of the matrix may occur. However, G-25 has been shown to withstand 0.1 M HCl for 1-2 hours and 0.02 M HCl for 6 months without any effect on its chromatographic properties.

Sephadex® resins are chemically resistant to 8 M urea. However, since such solutions would be very viscous, the flow rate would be much reduced in the presence of 8 M urea, and would lead to high back pressure. The beads are not able to withstand increased pressure to get a reasonable flow rate. In such cases, Sephacryl® resins should be used. Sephacryl® is more rigid and can withstand higher pressures. The Sephacryl® resin is also resistant to 8 M urea.

References

1. Porath, J., and Flodin, P., *Nature*, **183(4676)**, 1657-1659 (1959).
2. Stellwagen, E., *Meth. Enzymol.*, **182**, 317-328 (1990).
3. Ausubel, F.M. et al. (eds.), *Short Protocols in Molecular Biology*, 2nd edition. John Wiley & Sons (New York, NY), pp. 10-36 to 10-39 (1992).

MERCK

Notice

We provide information and advice to our customers on application technologies and regulatory matters to the best of our knowledge and ability, but without obligation or liability. Existing laws and regulations are to be observed in all cases by our customers. This also applies in respect to any rights of third parties. Our information and advice do not relieve our customers of their own responsibility for checking the suitability of our products for the envisaged purpose.

The information in this document is subject to change without notice and should not be construed as a commitment by the manufacturing or selling entity, or an affiliate. We assume no responsibility for any errors that may appear in this document.

Technical Assistance

Visit the tech service page at SigmaAldrich.com/techservice.

Standard Warranty

The applicable warranty for the products listed in this publication may be found at SigmaAldrich.com/terms.

Contact Information

For the location of the office nearest you, go to SigmaAldrich.com/offices.

The life science business of Merck operates as MilliporeSigma in the U.S. and Canada.

Merck and Sigma-Aldrich are trademarks of Merck KGaA, Darmstadt, Germany or its affiliates. All other trademarks are the property of their respective owners. Detailed information on trademarks is available via publicly accessible resources.

© 2021 Merck KGaA, Darmstadt, Germany and/or its affiliates. All Rights Reserved.

G2580pis Rev 09/21 MAM,SS,GCY

For research use only. Not for use in diagnostic procedures.

The Merck logo, consisting of the word "MERCK" in a bold, red, sans-serif font.