

Product Information

SIGMA QUALITY CONTROL TEST PROCEDURE Enzymatic Assay of **CHOLESTEROL OXIDASE**¹ (EC 1.1.3.6)

PRINCIPLE:

Cholesterol + O₂ $\xrightarrow{\text{Cholesterol Oxidase}}$ H₂O₂ + 4-Cholesten-3-one

H₂O₂ + o-Dianisidine (reduced) $\xrightarrow{\text{Peroxidase}}$ 2 H₂O + o-Dianisidine (oxidized)

CONDITIONS: T = 25°C, pH = 7.5, A_{500nm}, Light path = 1 cm

METHOD: Continuous Spectrophotometric Rate Determination

REAGENTS:

- A. 50 mM Potassium Phosphate Buffer, pH 7.5 at 25°C
(Prepare 100 ml in deionized water using Potassium Phosphate, Monobasic, Anhydrous, Sigma Prod. No. P-5379. Adjust to pH 7.5 at 25°C with 1 M KOH.)
- B. 1% (w/v) o-Dianisidine Solution (ODA)
(Dissolve the contents of a 50 mg vial of o-Dianisidine Dihydrochloride, Sigma Stock No. 510-50, in 5 ml of deionized water **or** prepare 5 ml in deionized water using o-Dianisidine Dihydrochloride, Sigma Prod. No. D-3252. **PREPARE FRESH.**)
- C. 0.5% (w/v) Cholesterol with 10% (v/v) Triton² X-100 Solution (Chol)
(Prepare by initially dissolving the Cholesterol, Sigma Prod. No. C-8503, in 10 ml of Triton² X-100. Heat until the solution clarifies. Then add 90 ml of deionized water. Stir well. Store the solution at 4°C.)³
- D. Peroxidase Enzyme Solution (POD)
(Immediately before use, prepare a solution containing 100 pyrogallol units/ml in deionized water using Peroxidase from Horseradish, Sigma Prod. No. P-8250.)
- E. Cholesterol Oxidase Enzyme Solution
(Immediately before use, prepare a solution containing 0.1 - 0.2 unit/ml of Cholesterol Oxidase in cold Reagent A.)

PROCEDURE:

Prepare a reaction cocktail by pipetting (in milliliters) the following reagents into a suitable container:

Reagent A (Buffer) 40.0
Reagent B (ODA) 0.50

Mix thoroughly by swirling and adjust to pH 7.5 at 25°C with 100 mM HCl or 100 mM KOH, if necessary. Then add Reagent A to a final volume of 50 ml. Mix by swirling thoroughly and oxygenate for approximately 10 minutes immediately before use.

Pipette (in milliliters) the following reagents into suitable cuvettes:

	<u>Test</u>	<u>Blank</u>
Reaction Cocktail	2.70	2.70
Reagent C (Chol)	0.10	0.10
Reagent D (POD)	0.10	0.10

Mix by inversion and equilibrate to 25°C. Monitor the A_{500nm} until constant, using a suitably thermostatted spectrophotometer. Then add:

Reagent A (Buffer)	-----	0.10
Reagent E (Enzyme Solution)	0.10	-----

Immediately mix by inversion and record the increase in A_{500nm} for approximately 5 minutes. Obtain the $\Delta A_{500nm}/\text{minute}$ using the maximum linear rate for both the Test and Blank.

CALCULATIONS:

$$\text{Units/ml enzyme} = \frac{(\Delta A_{500nm}/\text{min Test} - \Delta A_{500nm}/\text{min Blank})(3)(df)}{(7.5)(0.1)}$$

3 = Volume (in milliliters) of assay

df = Dilution factor

7.5 = Millimolar extinction coefficient of o-Dianisidine (Oxidized) at 500 nm

0.1 = Volume (in milliliters) of enzyme used

$$\text{Units/mg solid} = \frac{\text{units/ml enzyme}}{\text{mg solid/ml enzyme}}$$

$$\text{Units/mg protein} = \frac{\text{units/ml enzyme}}{\text{mg protein/enzyme}}$$

UNIT DEFINITION:

One unit will convert 1.0 μmole of cholesterol to 4-cholesten-3-one per minute at pH 7.5 at 25°C.

FINAL ASSAY CONCENTRATION:

In a 3.00 ml reaction mix, the final concentrations are 38 mM potassium phosphate, 0.009% (w/v) o-dianisidine, 0.017% (w/v) cholesterol, 0.33% (v/v) Triton X-100, 10 units peroxidase and 0.01 - 0.02 unit cholesterol oxidase.

REFERENCE:

Masurekar, P.S. and Goodhue, C.T. (1978) United States Patent No. 4,093,517, Eastman Kodak Company, Rochester, NY

NOTES:

1. This assay is not to be used to assay Cholesterol Oxidase from Schizophyllum commune, Sigma Prod. No. C-7274 and Cholesterol Oxidase from Brevibacterium sp., Sigma Prod. No. C-8153.
2. Triton is a registered trademark of Union Carbide Chemicals and Plastics Co., Inc.
3. The solution can still be used if it becomes cloudy upon addition of the water.
4. Peroxidase Unit Definition: One unit will form 1.0 mg purpurogallin from pyrogallol in 20 seconds at pH 6.0 at 20°C.
5. This assay is based on the cited reference.
6. Where Sigma Product or Stock numbers are specified, equivalent reagents may be substituted.

Sigma warrants that the above procedure information is currently utilized at Sigma and that Sigma products conform to the information in Sigma publications. Purchaser must determine the suitability of the information and products for its particular use. Upon purchase of Sigma products, see reverse side of invoice or packing slip for additional terms and conditions of sale.