

Product Information

4-Nitrophenyl phosphate disodium salt hexahydrate

Tablet (15 mg substrate per tablet)

N2640

Product Description

CAS Registry Number: 333338-18-4

Synonyms: pNPP, *p*-Nitrophenyl Phosphate, Phosphatase substrateMolecular Formula: $C_6H_4NNa_2O_6P \cdot 6 H_2O$

Formula Weight: 371.14

4-Nitrophenyl Phosphate (pNPP) is the substrate of choice for use with alkaline phosphatase conjugates in Enzyme Linked Immunosorbant Assay (ELISA) procedures, due to its high sensitivity.^{1,2} ELISA applications that use pNPP may be read in timed assays or stopped with alkaline solutions for delayed readings.³ This substrate produces a soluble end product that is yellow in color and can be read spectrophotometrically at 405 nm. The pNPP reaction may be stopped with the addition of 3 M NaOH solution and read at 405 nm.

The N2640 tablets are supplied as 50 tablets (50TAB) or 100 tablets (100TAB) per box, individually foil wrapped for ease of use, storage, and safety. Several publications,⁴⁻⁶ theses⁷ and dissertations⁸⁻¹² have cited use of N2640 in their research protocols.

Precautions and Disclaimer

This product is for R&D use only. Not for drug, household, or other uses. Please consult the Safety Data Sheet for information regarding hazards and safe handling practices.

Storage/Stability

The tablets should be stored at -20 °C.

Preparation Instructions

Tablets should be warmed to room temperature before use. Solutions should be freshly prepared.

Dissolve 1 tablet in either Glycine Buffer or Diethanolamine Buffer to the desired concentration. (Typically a pNPP concentration of 1 mg/mL is used.)

Glycine Buffer

To prepare 0.1 M glycine buffer, pH 10.4, with 1 mM $MgCl_2$ and 1 mM $ZnCl_2$:

- Add 7.51 g of glycine, 203 mg of $MgCl_2$, and 136 mg of $ZnCl_2$ to ~980 mL of water. Mix.
- Adjust pH to 10.4 with 19 M NaOH.
- Bring the volume to 1 L with water.

Diethanolamine Buffer

To prepare 1 M diethanolamine buffer, pH 9.8, with 0.5 mM $MgCl_2$:

- Add 97 mL of diethanolamine and 100 mg of $MgCl_2$ to 800 mL of water.
- Adjust pH to 9.8 with 10 M HCl.
- Bring the volume to 1 L with water.

Procedure

General ELISA procedure with alkaline phosphatase conjugates

- Add 200 μ L of substrate solution (typically 1 mg/mL) per well.
- Incubate the plate in the dark for 30 minutes at room temperature.
- The absorbance can be read at 405 nm on a multiwell plate reader.
- The reaction may be stopped by adding 50 μ L of 3 M NaOH per 200 μ L of reaction mixture.

Troubleshooting

If the background is too high:

1. Use a blocking step prior to the application of the primary antibody. Normal Serum (5% v/v) from the same species as the host of the second antibody generally produces the best results.
2. Additional blocking agents for an ELISA are:
 - 0.05% TWEEN® 20 in 50 mM TBS, pH 8.0.
 - 1% BSA containing 0.05% TWEEN® 20 in 50 mM TBS, pH 8.0.

- 3% nonfat-dried milk in 0.01 M TBS (Cat. No. P2194). **Do not use milk as a blocking agent when using avidin-biotin systems.**
3. Use 0.05% TWEEN® 20 in all washing and antibody diluent buffers.
 4. Run control wells without the primary antibody to check for non-specific reactivity of the secondary antibody/alkaline phosphatase conjugate.
 5. Adjust the titer of the primary antibody and/or the alkaline phosphatase conjugate to determine the optimal working dilutions.

If no color develops or color is too faint:

1. Adjust the concentration of the primary antibody.
2. Adjust the concentration of the secondary antibody/alkaline phosphatase conjugate.
3. Determine if the enzyme conjugate is active by mixing a small sample of substrate and conjugate together in a test tube.
4. Increase the substrate incubation time or temperature.
5. Adjust the concentration of the coating antigen.
6. Consider using an amplifying system such as avidin-biotin.

References

1. Voller, A. et al., *Bull. World Health Organ.*, **53(1)**, 55-65 (1976).
2. Engvall, E., *Methods Enzymol.*, **70(A)**, 419-439 (1980).
3. Voller, A., and Bidwell, D., "Enzyme-linked immunosorbent assay", in *Manual of Clinical Laboratory Immunology*, 3rd ed. (Rose, N.R. et al., eds.). American Society for Microbiology (Washington, D.C.), pp. 99-109 (1986).
4. Malla, R. et al., "Immuno-Technology for the Localization of Acid Phosphatase Using Native Gel Bands in *Piriformospora indica* and Other Soil Microorganisms", in *Advanced Techniques in Soil Microbiology* (Varma, A., and Oelmüller, R., eds.). Springer-Verlag (Berlin), pp. 213-235 (2007).
5. Sample, C.J. et al., *Peptides*, **48**, 96-105 (2013).
6. Surman, S.L. et al., *Clin. Vaccine Immunol.*, **21(4)**, 598-601 (2014).
7. Barefoot, Brice Edward, "Increasing the Immunogenicity of Vaccinia Virus Vectors by Insertion of Cowpox Virus A-Type Inclusions". North Carolina State University, M.S. thesis, p. 36 (2012).

8. Furrie, Elizabeth, "Intestinal modification of a protein antigen and its effect on oral tolerance induction". University of London, Ph.D. dissertation, p. 44 (1993).
9. Zakzuk, Josefina, "Immunological and Biological Characterization of *Blomia tropicalis* Allergens". University of Cartagena, Ph.D. dissertation, p. 172 (2012).
10. Ausländer, David, "Engineering of synthetic biology-based gene networks in mammalian cells for biomedical applications". ETH Zürich, Dr. sc. dissertation, p. 161 (2014).
11. Lahl, Annika Olivia Regina, "Der Einfluss von Vitamin D auf die spezifische humorale Immunantwort" ("The influence of vitamin D on the specific humoral immune response"). Charité – Universitätsmedizin Berlin, Dr. med. dissertation, p. 77 (2015).
12. Tastanova, Aizhan, "Synthetic Mammalian Gene Regulation Systems for Biomedical Applications". ETH Zürich, Dr. sc. dissertation, p. 39 (2016).

Notice

We provide information and advice to our customers on application technologies and regulatory matters to the best of our knowledge and ability, but without obligation or liability. Existing laws and regulations are to be observed in all cases by our customers. This also applies in respect to any rights of third parties. Our information and advice do not relieve our customers of their own responsibility for checking the suitability of our products for the envisaged purpose.

The information in this document is subject to change without notice and should not be construed as a commitment by the manufacturing or selling entity, or an affiliate. We assume no responsibility for any errors that may appear in this document.

Technical Assistance

Visit the tech service page at [SigmaAldrich.com/techservice](https://www.sigmaaldrich.com/techservice).

Terms and Conditions of Sale

Warranty, use restrictions, and other conditions of sale may be found at [SigmaAldrich.com/terms](https://www.sigmaaldrich.com/terms).

Contact Information

For the location of the office nearest you, go to [SigmaAldrich.com/offices](https://www.sigmaaldrich.com/offices).

The life science business of Merck operates as MilliporeSigma in the U.S. and Canada.

Merck and Sigma-Aldrich are trademarks of Merck KGaA, Darmstadt, Germany or its affiliates. All other trademarks are the property of their respective owners. Detailed information on trademarks is available via publicly accessible resources.

© 2013-2023 Merck KGaA, Darmstadt, Germany and/or its affiliates. All Rights Reserved.
N2640pis Rev 02/23 RBG,CMH,MAM,GCY

